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JM109(DE3) Chemically Competent Cell

Cat.NO. ZC1230

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Sample Induction Protocol (for reference only)

1. Inoculate a single colony from a freshly streaked plate into 3ml of LB medium containing the appropriate antibiotic for the plasmid and host strain.
2. Incubate with shaking at 200rpm at 37°C overnight.
3. Inoculate 50ml of LB medium containing the appropriate antibiotic with 0.5ml of the overnight culture prepared in step 2 (use the 500ml triangular flask as the container would be better).
4. Incubate with shaking at 150rpm at 37°C until the OD 600 reaches 0.5-0.8. (0.6 recommended; about 2.5h).
5. (Optional) Pipet 1ml of the cultures into clean microcentrifuge tubes and place the tubes on ice until needed for gel analysis or storage at -20°C. These will serve as the non-induced control samples.
6. Add IPTG to a final concentration of 1mM. Optimal time for induction of the target protein may vary from 2-16 hours, depending on the protein.
7. Incubate with shaking at 120rpm at 37°C for 2-4 hours. To determine the optimal time for induction of the target protein, it is recommended that a time course experiment be performed varying the induction from 2-16 hours.
8. Place the culture on ice for 10 minutes. Harvest cells by centrifugation at 5,000 × g for 10 minutes at 4°C.
9. Remove the supernatant and store the cell pellet at -20°C (storage at lower temperatures is also acceptable).

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